Operation Manual V1.1





This Product is for Research Use Only



Thank you for purchasing the Smart-Q[™] 800 Fluorometer. This manual details the instrument's features, specifications, as well as complete operating instructions; please read it carefully prior to operation. Keep this user manual for later use.

Important:

Please keep the box and packaging material for this instrument. If service is required, the box will be needed to ship the instrument to our Service Department.

Initial Inspection

Please inspect the instrument as well as all included accessories when you first open the packaging. If you find anything damaged or missing, please contact Benchmark Scientific or your local distributor immediately.

BENCHMARK SCIENTIFIC / ACCURIS INSTRUMENTS

PO Box 709 Edison, NJ 08818 USA Phone: 908-769-5555 Website: www.benchmarkscientific.com/ www.accuris-usa.com Email: info@benchmarkscientific.com / info@accuris-usa.com

Safety Warning and Guidelines

1. Important information for safe use

Users should have a clear understanding of how to use this instrument before operation, please read this manual carefully prior to operation.



NOTICE: Please review this manual in its entirety before operating the Smart-Q[™] 800.



For Research Use Only. This instrument is not a medical device and is not intended to be used for clinical testing.

2. Operation and Maintenance

The operation and maintenance of the instrument should comply with the basic guidelines and warnings below. Incorrect operation or maintenance will have detrimental effects on the life, performance, and safety features of the instrument.



This instrument is for indoor use only and conforms to I class B type common equipment of standard GB9706.1.



CAUTION: Biological Contamination

All samples used with this instrument, including quality control samples, calibration samples, and components should be considered infectious and treated with caution. Always wear gloves and appropriate safety gear for protection.



The instrument should not be disassembled by the end user unless under the guidance of Benchmark's Service Department.



The instrument should only be installed in an environment with standard room temperature, low dust, low humidity, and away from direct sunlight, electromagnetic interference, and heat sources.



Always power off the instrument when you are finished using it. Unplug the power cord and cover the instrument with a cloth or plastic sheet to prevent excessive dust from entering the housing.



Unplug the instrument immediately if any of the following occur, and contact Benchmark Scientific's Service Department:

- Liquid enters the Instrument.
- The instrument is dropped or the outer shell damaged.
- The functions of the instrument have obviously changed.

3. Maintenance

The instrument housing should be cleaned regularly using a soft, damp cloth.

To clean the touchscreen, disconnect the power cable, and clean the touchscreen with a soft cloth lightly moistened with LCD cleansing detergent. Take care not to use excessive force as this can damage the touchscreen. Do not use abrasive cleaning solutions or material to prevent the touchscreen from getting scratched.

To disinfect the instrument, disconnect the power cable and use a soft cloth lightly moistened with 70% ethanol, 70% isopropanol, or 10% bleach (0.6% sodium hypochlorite). Take care so that the cleaning solution does not enter the power inlet, the sample port, or the USB drive ports.

4. Transportation and Storage Requirements

Ambient temperature: 10°C ~ 40°C Relative humidity: ≤ 70% Place in a well-ventilated room, away from corrosive gas.

Table of Contents

Chapte	1 Introduction
Chapte	2 Specifications
1.	Required Installation Environment4
Env	ironmental Temperature: 10°C~40°C4
2.	Specifications4
Chapte	3 Instrument Overview
Chapte	4 Installation
4.1	Initial Check6
4.2	Installation Conditions6
4.3	Operating Conditions6
Chapte	5 Software Introduction
5.1	Instrument Startup8
5.2	Main Menu Interface8
5.3	dsDNA Interface9
5.3	1 dsDNA Interface10
5.3	2 Calibration Curve Interface
5.3	3 Sample Measurements11
5.3	4 Molarity Calculator
5.1	1 Normalization Calculator13
5.2	Fluorescence Interface14
5.3	Working Solution Calculation Settings15
5.4	Standard Curve Interface16
5.4	1 New Standard Curve16
5.4	2 Curve Management
5.5	Reports Interface18
5.6	User Management Interface19

Ch	apter 6 T	roubleshooting	22
	5.7.3	Maintenance Interface	21
	5.7.2	Upgrade Interface	21
	5.7.1	System Settings Interface	20
	5.7	Settings Interface	20

Chapter 1 Introduction

The Smart-Q[™] 800 Fluorometer is used for the precise quantitation of nucleic acids and proteins in microliter volumes via fluorescent detection. Under conditions of low concentrations of molecules, concentration values are linear to fluorescence intensity. Thus, unknown samples can be quantitatively analyzed. Reagent kits with fluorescent dye, buffer, and standards are available from Accuris or your local distributor.

Key Features and Benefits:

- > 7-inch color LCD touch screen display, easy to operate
- Measure up to 8 samples at a time with 8-strip tubes (0.2mL)
- Fast and accurate detection of DNA, RNA and proteins (< 6 sec)</p>
- Low sample consumption, only 1-20µL of sample is needed for accurate quantification
- Data storage for up to 10,000 sample records
- Sample data can be printed or exported via USB drive
- Customized sample standard curve
- Reagent calculator to assist with volume determinations
- Auto sleep function for power-saving

Chapter 2 Specifications

1. Required Installation Environment

Environmental Temperature: 10°C~40°C Relative Humidity: ≤ 70% Input Voltage: 24 VDC 2A

2. Specifications

Model	Smart-Q [™] Fluorometer
Display	7.0-inch, color touch screen
Dynamic Range	4 orders of magnitude
Light Course	Blue LED
Light Source	Red LED
Evolution Woyalangth	Blue: 470 ± 15nm
Excitation Wavelength	Red: 625 ± 20nm
Emission Wayslongth	Green: 515 – 540nm
Emission Wavelength	Red: 670 – 725nm
Stability	≤ 1.5%
Detector	Photodiodes (300-1100nm)
Measurement Speed	≤ 6 seconds
Calibration type	2-3-point standard
Consumable type	8-strip optically clear PCR tubes (0.2mL)
Input Voltage	24 VDC 2A
Dimensions (WxDxH)	16.1 x 28.7 x 7.6 cm / 6.3 x 11.3 x 3.0 in.
Weight	2.0kg / 4.4lbs

Chapter 3 Instrument Overview

This chapter provides an overview of the physical characteristics of the Smart-Q 800. When using the instrument for the first time, please read this chapter carefully to familiarize yourself with all the parts.

Structure 1:

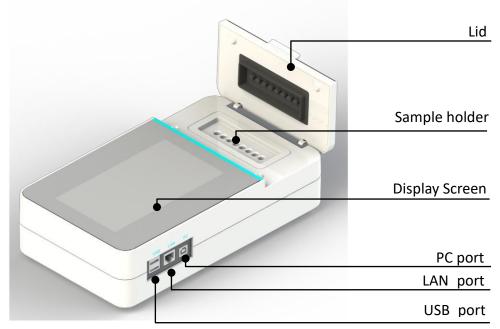


Fig 3.1.01 Front/Side View

Structure 2:

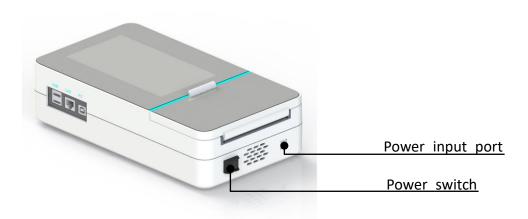


Fig 3.1.02 Rear View

Chapter 4 Installation

4.1 Initial Check

Each Smart-Q 800 has been factory-inspected prior to shipping. Please refer to the included packing list to ensure that all items are included. Contact Accuris or your local distributor if there are any signs of damage to the packaging upon receipt.

Carefully check the instrument's appearance and housing for cracks, damage, or deformation.

4.2 Installation Conditions

Install the instrument on a clean, dry, and level surface in a well-ventilated room away from corrosive agents

Installation Steps

- 1. Carefully remove the instrument from its packaging and place it onto a stable worktable. Remove all plastic and tape.
- Unpack the 24V 2A power adapter and plug it into the socket on the back of the instrument. Then connect another side to an AC100~240V power outlet.
- 3. Turn on the power. The instrument will start to self-check after powering on.

4.3 Operating Conditions

The Smart-Q uses an 8-strip (0.2ml) fluorescent PCR tube to test samples.

Sample test steps:

1 Insert the 0.2mL tube with sample or standards into sample holder.

2 Close the lid; Click the measurement button on the interface to perform a measurement.

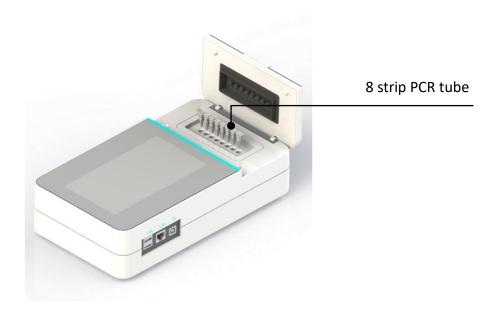


Fig 4.3.01 Instrument Operation Diagram

Chapter 5 Software Introduction

5.1 Instrument Startup

Power the instrument on. Upon start-up, the instrument will perform a self-check (**Fig. 5.1.01**). After the self-check, the Login Interface will be displayed (**Fig. 5.1.02**).



Fig 5.1.01 Self-check



Fig 5.1.02 Login Interface

5.2 Main Menu Interface

After entering the user-specific password (**Admin password is set by default to** "0000"), the Main Menu Interface will be displayed (**Fig. 5.2.01**).



Fig 5.2.01 Main Menu Interface

Name	Function
	Select to manage account settings or to switch between users.
\gg	Select to view the working solution calculator
diDNA	Select to enter the dsDNA Interface
ji) ¹ RIVA	Select to enter the RNA Interface
Oligo	Select to enter the Oligo Interface
Protein	Select to enter the Protein Interface
Fluorescence	Select to enter the Fluorescence Interface
Curve	Select to enter the Standard Curve Interface
Report	Select to view a list of the most recent protocols created/run
Settings	Select to enter the settings interface

Table 5-1 Main Menu Interface Operations

5.3 dsDNA Interface

Note: As the interface of dsDNA, RNA, Protein, Oligo are identical, Only the dsDNA Interface is described in detail in this manual.

From the Main Menu Interface, click to enter the dsDNA Interface. Three test types are available for selection: dsDNA: High Sensitivity, dsDNA: Broad Range, & 1x dsDNA (**Fig. 5.3.01**).

×	<	dsDNA: High Sensitivity
Admin 03/18/2022 09:51:45	DNA-1243	16 🖉
	Curve name	dsDNA01 🔻 🔟
and the second s	Unit	ng/µL ▼
Test type		
dsDNA: High Sensitivity	S1 S	2 53 54 55 56 57 58
dsDNA: Broad range		
🔿 1x dsDNA	Volume(j	il) 10
Cancel OK		
Curve Report Settings		$\langle \rangle$
		•• →
Fig 5.3.01 dsDNA test type	Fig 5	.3.02 dsDNA test

5.3.1 dsDNA Interface

Select the test type and click "OK" to enter next interface.

The following takes *dsDNA: High Sensitivity* as an example.

The specific functions are as follows:

ng/µL

Curve name dsdnaCurve02 : Select an available standard curve.

Sample name S1 : Input the sample name.

 \mathbf{I} : Select the appropriate units: ng/µL, ng/mL, µg/µL, µg/mL,

mg/mL.

Unit

Click to enter the Calibration Curve Interface (see Chapter 5.3.2).

: Click to print all current protocol test data.



: Click to perform a sample measurement.



: Click to modify the protocol name.

Volume(µl)

10

: Set the volume of the original sample, range 1~20µL.

Original Conc.: Calculated concentration of the original sample.

RFU (Relative Fluorescence Units): Calculated fluorescence value of the original sample.

5.3.2 Calibration Curve Interface

Prior to sample measurements, a calibration curve must be created using the standards that are typically included with a quantification kit. Only 2 standards are required to form the calibration curve.

Enter the dsDNA testing interface. Click the button and select "Calibration" to create the calibration curve (**Fig. 5.3.03**).

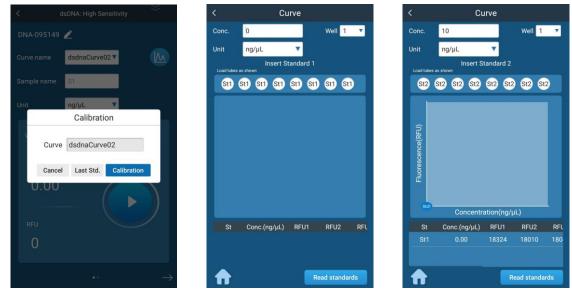


Fig 5.3.03 Curve Cal.

Fig 5.3.04 First Cal. point

Fig 5.3.05 Second Cal. point


For each assay, you have the choice to run new standards, or to use values from the previous calibration to calibrate the assay. If a calibration has been carried out recently, the last set data can be selected; If recalibration is required, select the button to perform the calibration. Up to 8 of each standard can be used for calibration. To select certain wells for calibration (i.e., Well 1 and 2 only), select set to enable the specified well. The "St1" icon will turn blue in color once enabled. **Note:** Samples can only be measured in wells for which calibration has been performed; For example: if calibration was performed on wells 1 and 2, then samples can only be measured in wells 1 and 2.

Enter the calibration interface, enter the known concentration value for each standard, and press "Read Standards" to perform measurements of the standards. After both standards have been read, select "Save" to save the calibration data.

5.3.3 Sample Measurements

After generating the calibration curve, input the original volume of the sample that was added to the tube and click is to perform a measurement. After measurement, the original concentration and RFU of the sample will be displayed. Swipe right to view

data from multiple samples within the same run. Click the name of the sample to modify the sample's name.

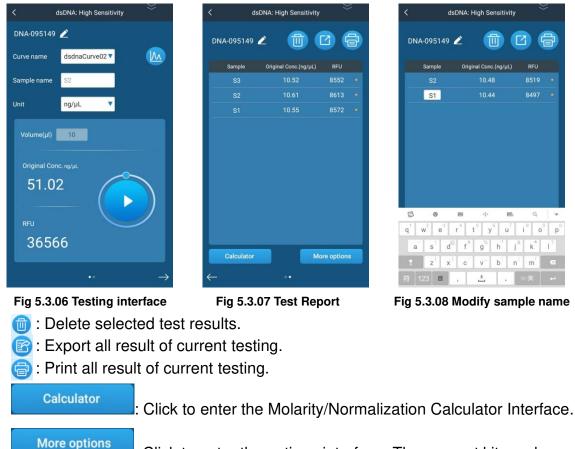

: Click to enter the options interface. The reagent kit number and sample ID can be adjusted in this interface.

5.3.4 **Molarity Calculator**

Click

Calculator

and select the "Molarity" tab to enter the Molarity interface.



Fig 5.3.09 Molarity/Normalization Calculator Interface



Select the conversion units: $ng/\mu L$, ng/mL, $\mu g/\mu L$, $\mu g/mL$, mg/mL.

After selecting the conversion units, select the units: M, mM, µM, nM, pM.



Input the molecular weight; the default value for dsDNA is set to 660g/mol.



Input the DNA length, when "Auto-populate DNA length" in enabled, all lengths will be automatically filled in according to the length of the first sample.

Note: The Molarity/Normalization Calculator Interface can only be selected one sample measurements have been performed.

5.1.1 Normalization Calculator

Select the Normalization tab to enter the Normalization calculator interface (**Fig. 5.3.10**). In this interface, choose the type of Normalization (for example: molarity, concentration, and/or weight), and set the final sample molarity and volume. Click "Calculate" to enter the results interface (**Fig. 5.3.11**).



Fig 5.3.10 Normalization Interface



Fig 5.3.11 Results Interface

The Filling List tab shows the required sample and buffer volume to be added to achieve the desired final sample concentration within the final sample volume. On the Conc. Ratio interface, if a secondary dilution is required, the dilution ratio and the concentration after dilution will be displayed (**Fig. 5.3.12**). The Initial Conc. interface shows the concentration of the original sample (**Fig. 5.3.13**). Click "Export" to export current data to an inserted USB drive.

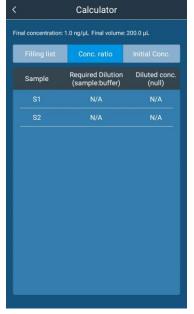


Fig 5.3.12 Secondary Dilution

inal concentration: 1.0 ng/µL Final volume: 200.0 µL Filling list Conc. ratio Initial Conc Sample Concentration(ng/µL) S1 51.02 S2 51.00	<		Calculator	
Sample Concentration(ng/µL)	Fin	al concentration	n: 1.0 ng/µL. Final volum	ie: 200.0 µL
S1 51.02		Filling list	Conc. ratio	Initial Conc
		Sample	Concentration(ng/	μL)
S2 51.00			51.02	
			51.00	

Fig 5.3.13 Initial Concentration of Sample

5.2 Fluorescence Interface

From the Main Menu Interface, select it to enter the Fluorescence interface. Three test types are available for selection: Blue 470nm, Red 635nm, & Blue & Red (**Fig. 5.4.01**).

Note: The Fluorescence interface does not require the generation of a calibration curve.

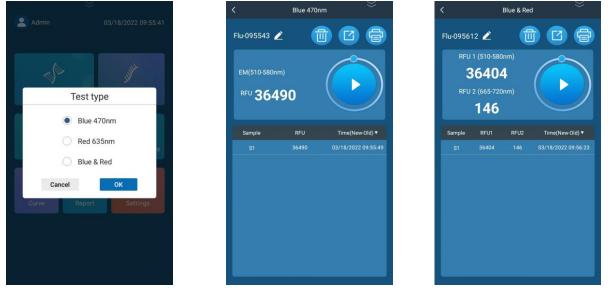


Fig 5.4.01 Fluorescence Test Type

Fig 5.4.02 Blue 470nm Channel

Fig 5.4.03 Blue & Red Channels

Test Types: Blue 470nm and Red 635nm are single excitation wavelength tests, whereas Blue & Red is a dual-wavelength test (**Fig 5.4.03**).

The Blue 470nm interface is identical to the Red 635nm interface. Click to begin a sample measurement. After measurement, the RFU of the sample will be displayed (**Fig. 5.4.02**). Data can be deleted, exported, or saved via the icons found at the upper right corner of the display.

5.3 Working Solution Calculation Settings

Click and drag down the button to view the Working Solution Calculator (**Fig. 5.5.01**). This interface allows for the calculation of the total working solution volume required based on dye-to-buffer ratios. Input the number of samples and standards required and select "Done". The calculated volumes for "Dye", "Buffer", & "Total volume" will be displayed based on the set numbers of samples and standards. Select "Include Overage" to account for 1 extra tube in the calculation.

Detection Range Calculator: From the Working Solution Calculator interface, swipe right to enter the Detection Range Calculator (**Fig. 5.5.02**). Here, the assay type, units, and original sample volume can be adjusted, and a detection range will be displayed after clicking "Done".



Fig 5.5.01 Working Solution Calculator

Fig 5.5.02 Detection Range Calculator

to enter the Standard Curve Interface (Fig.

5.4 Standard Curve Interface

From the Main Menu Interface, click **5.6.01**).

All
Type

Name *
Type

Name *
Type

Mame *
Type

MacOurve33
RNA ER

MacOurve34
RNA ER

MacOurve24
Micro RNA

Users
MacOurve22


RNA ER
Users

macOurve21
RNA HS

Users
MacOurve21

MacOurve21
RNA HS

Fig 5.6.01 Standard Curve Interface

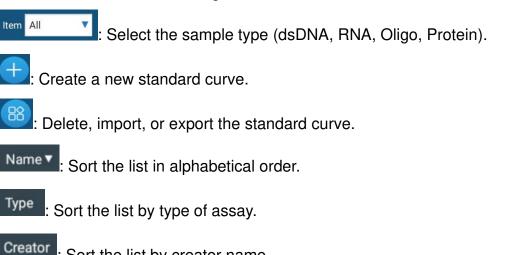

: Sort the list by creator name.

5.4.1 New Standard Curve


Click the \bigcirc icon to create a new standard curve. The New Curve Selection Box will be displayed, with adjustable values for sample type, assay type, and curve name (**Fig. 5.6.02**). Once the values have been adjusted/set, click "OK" to enter the New Standard Curve Interface (**Fig. 5.6.03**).

<		Curve		
Item All		2		
		New Curve	9	
	dsDNA		~	
	dsDNA	High Sensitiv	ty 🗸	
	DNA09	5926		
	Cance		ок	
maCur	ve22	RNABR	Users	
			/2 >	1

Fig 5.6.02 New Curve Selection Box



Fig 5.6.03 New Standard Curve Interface

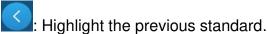

Highlight the next standard, or to add an additional standard.

- $^{
 u}$: View Parameter Settings (EX / EM wavelengths + concentration units).
 - $\stackrel{\frown}{=}$: Plot the obtained sample data points to a curve.
- Uzeration: Delete RFU1/RFU2/RFU3 data or to delete the standard itself.
 - : Perform standard sample measurement.

Note: At least 5 standards are required to create the standard curve. For each standard, the average RFU value taken from 3 readings is used for curve generation.

5.4.2 Curve Management

From the Standard Curve Interface Click 88 to enter Curve Management Interface

(Fig.	5.6.04)
-------	-----------------

圃

Г7

Cancel	Curve	
Item All	• 💼	
Name ▼	Туре	Creator
maCurve34		Tester
maCurve33	RNA ER	Tester
maCurve32	RNA BR	Tester
maCurve31	RNA HS	Tester
maCurve24		
maCurve23	RNA ER	
maCurve22	RNA BR	
maCurve21		
maCurve14		
Total: 27	< 1/2	2 > 1

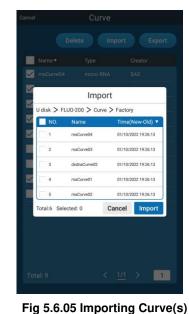


Fig 5.6.04 Curve Management Interface

- Delete the selected curve.
- Export the selected curve(s) to a USB drive (Fig 5.6.04).
- : Import curve(s) from a USB drive to the instrument (Fig. 5.6.05).

5.5 Reports Interface

In the main interface, click is to enter the Reports Interface (**Fig. 5.7.01**). Input keywords in the search box to quickly find the required report (**Fig 5.7.02**).

	Report	
2 Keyword		8
Name	Curve	Time(New-Old)
		03/18/2022 09:56:12
		03/18/2022 09:55:57
		03/18/2022 09:55:43
DNA-095149	dsdnaCurve02	03/18/2022 09:51:49
Flu-093740		03/18/2022 09:37:40
Flu-093724		03/18/2022 09:37:24
		03/18/2022 09:37:13
		03/18/2022 09:36:59
Flu-093648		03/18/2022 09:36:48
Total: 66		5 > 1

Fig 5.7.01 Reports Interface

<	Report	
Q dna	0	88
Name	Curve	Time(New-Old) ▼
DNA-095149	dsdnaCurve02	03/18/2022 09:51:49
DNA-093119	dsdnaCurve02	03/18/2022 09:31:19
DNA-092633	dsdnaCurve02	03/18/2022 09:26:33
DNA-092043	dsdnaCurve02	03/18/2022 09:20:43
	dsdnaCurve02	03/17/2022 15:20:52
	dsdnaCurve02	03/17/2022 14:15:29
DNA-133845	dsdnaCurve01	03/17/2022
8 0	四 ·I·	₩ Q -
$\mathbf{q}^{\dagger} \mathbf{w}^2 \mathbf{e}^3$	r ⁴ t ⁵ y ⁶	u ⁷ i ⁸ o ⁹ p ⁰
a s d	f [#] g [%] h	$j^{\delta} k^{\star} l^{?}$
1 z (x)	c v b	n' m' 🗠
符 123 回	, 🛓	. 中/英 ←

Fig 5.7.02 Quick Search

Click on a report to view its saved data (**Fig. 5.7.03**). Click on the sample from the row to view the detailed sample report (**Fig. 5.7.04**).

<	DNA-090433	
 Test type 	1x dsDNA	
Volume	10.0µL	
Operator	Admin	
Excitation	Blue	
Time	03/15/2022 09:04:33	
Standard curve	dsdnaCurve11	
Creator	Admin	
Standard 1	Conc.3.00ng/µL RFU 3.3	
Standard 2	Conc.9.00ng/µL RFU 9.9	
Calibration time	03/15/2022 10:04:24	
Sample	Original Conc.(ng/µL)	RFU
S1		
S2		
S3		
S4		
S5		
S6		
S8		

Calculator	Delete	Export

Fig 5.7.03 Saved Report Data

S1 Info.	
Original concentration	:1.10ng/µL
Sample RFU value	:1.3
Molarity calculation I	nfo.
Desired unit	:ng/µL->nM
Molecular weight	:660
Length(bp)	:6000
Molarity	:3.00nM
Normalization calcula	ation Info.
Final concentration	:1.00ng/mL
Final volume	:200µL
Add sample	:1.09µL
Add buffer	:198.91µL
Required Dilution	:1:5
Diluted concentration	:183.33ng/mL
Concentration	:1100.00ng/mL

Fig 5.7.04 Detailed Sample Report

5.6 User Management Interface

From the Main Menu Interface, select the Sicon from the upper left corner of the display. The following icons will be displayed:

(User Settings: Administrators can manage user accounts by adding, deleting, or modifying the user name / password. Non-Administrators can only change their own passwords in this interface.

=) : Switch User: Return to the Login Interface to switch to a different user.

5.7 Settings Interface

Click

in the main interface to enter the Settings Interface (Fig. 5.9.01).



Fig 5.9.01 Settings Interface



5.7.1 System Settings Interface

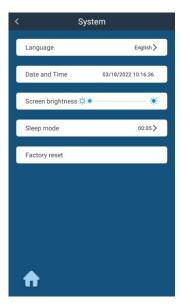


Fig 5.9.02 System Settings

to enter system settings (Fig. 5.9.02):

Language: Select English, or Chinese.

Click

Date and Time: Set the system date and time.

Screen Brightness: Adjust the screen brightness.

Sleep Mode: Turn the automatic sleep mode on or off and adjust time required to initiate Sleep Mode.

5.7.2 Upgrade Interface

From the Main Menu Interface, click \checkmark to enter the Upgrade Interface (**Fig 5.9.03**). The versions of the software and firmware will be displayed. To make updates, insert a USB drive with the relevant update files and select the ">" next to either the software or firmware version number.



Fig 5.9.03 Upgrade interface

5.7.3 Maintenance Interface

From the Main Menu Interface, click and input the password to enter the Maintenance Interface. This is only available for authorized service procedures. Contact Accuris Instruments or your local distributor for details.

Chapter 6 Troubleshooting

No.	Issue	Analysis	Solutions
1	No display after turning on power	Power not reaching the instrument	Check the power supply; Check whether the plug is loose.
2	Light source not visible	Light source failure; Module poor contact	Check the power; Check whether the module is loose, if loose Contact Accuris Instrument or your local distributor
3	Measurement not stable	PCR strip isn't in place	Check the PCR strip and press it in place
4	Measurement not accurate	Standard curve is not accurate; Instrument detection is drifting	Check standard solution and reset standard curve; Calibrate standard curve

Quantification Kits and Detection Tubes



NS1020-HS-100 High Sensitivity DNA Quantification Kit, 100 assays	
NS1020-BR-100	Broad Range dsDNA Quantification Kit, 100 assays
NS1020-T	0.5mL Fluorescence Detection Tubes, pack of 250