



Electrophoresis Power Supply



- Touch screen control Simple programming and operation
- Bright color display
 Shows all parameters simultaneously
- Dual orientation design Fits on a crowded bench or shelf
- High performance specifications

Max Amperage:

300 V 400 mA

Max Wattage:

Max Voltage:

100 Watts





Touch Screen

Rotating Display

Compatible With:







Horizontal **Tanks**



CE

Blotting Systems



Exclusive "Dual Orientation" feature (saves lab space)



Powerful and Capable

The Accuris myVolt™Touch is a work horse electrophoresis power supply, ideal for efficient separation of nucleic acids and proteins as well as basic western blotting applications. Four pairs of output jacks are wired in parallel for connecting multiple electrophoresis tanks. A color touch screen allows quick setting and monitoring of all run conditions, including run time, voltage, and current. Users can select "Constant Current" or "Constant Voltage" during an electrophoresis run, keeping the chosen parameter fixed as resistance changes. The dual orientation design with rotating display allows the myVolt Touch to be installed in the most convenient location, on a shelf or bench, saving valuable space.

Touch Screen Control Panel:

The myVolt Touch features a logical user interface. Setting up an electrophoresis run couldn't be easier; simply touch the parameter to select it, press the "up" and "down" arrows to adjust the value, and then touch "Start". All parameters are always displayed, with no need to toggle between Time, Voltage or Amperage. In addition, the run status is always displayed.

Time:	01:00	:00
Voltag	je: 10	0 v
mAmı	: 400	mA
Sta	atus: ST	OPPED
Start		Constant

Specifications		
Output Settings:	Constant Voltage or Constant Current	
Voltage Output:	10 to 300V (set in 1V increments)	
Current Output:	4-400mA (set in 1mA increments)	
Max Wattage:	100W	
Output Terminals:	4 pairs, in parallel	
Timer:	1 min 20 hrs, or Continuous	
Safety:	No load detection, overload detection with power limiting function	
Display:	2x2 in. color touchscreen	
Weight:	2.0 lbs./0.9 kg.	
Dimensions:	8.5x6.3x4.0in. (21.5x15.5x9.75cm)	
Electrical:	100V to 230V, 50-60Hz (Universal Input)	

Ordering Information:

E2301* myVolt Touch Power Supply, 115V

*115V with US plug. For 230V, please add (-E)

Also available from Accuris:





EM: info@accuris-usa.com