

		<u>DNA Stain Type</u>	
		Ethidium Bromide	SmartGlow™ SYBR® Green
<u>Excitation Source</u>	SmartDoc blue illumination base	Not recommended	E5001-ORANGE Orange PMMA Filter
	Blue light transilluminator	E5001-ORANGE Orange PMMA Filter	E5001-ORANGE Orange PMMA Filter
	UV transilluminator	E5001-590 590nm Filter	E5001-535 535nm Filter

For optimum image quality:

- Remove the phone's case
- Line up the camera lens with the photo filter

Place the photo filter into the recess on the top platform

Remove the extender for close ups or small gels



E5001-ORANGE Orange PMMA Filter (use with blue light)



E5001-535 535nm Filter (use with UV)



E5001-590 535nm Filter (use with UV)